

Instruction Manual

DEMI-02

DEMI-05

DEMI-10

Demineralizer Bottle



Dear Customer,

Thank you for choosing a Hanna Product.

Please read the instruction sheet carefully before using the chemical kit. It will provide you with the necessary information for correct use of the kit. If you need additional information, do not hesitate to e-mail us at tech@hannainst.com.

Remove the chemical kit from the packing material and examine it carefully to make sure that no damage has occurred during shipping. If there is any noticeable damage, notify your Dealer or the nearest Hanna office immediately.

Each kit is supplied with:

- Demineralizer Bottle with filter cap;
- 1 calibrated plastic beaker (100 ml);
- Instruction Manual

Note: Any damaged or defective item must be returned in its original packing materials.

SPECIFICATIONS

Code Kit	for preparation of.....DI water
DEMI-02	2 liters
DEMI-05	5 liters
DEMI-10	10 liters

Note: The exact amount of water that can be demineralized with a single Demineralizer Bottle varies depending on the hardness level of the tap water.

HEALTH AND SAFETY

The chemicals contained in this kit may be hazardous if improperly handled. Read the relevant Health and Safety Data Sheets before using the kit.

GENERAL DESCRIPTION

The Demineralizer Bottle contains a mixed bed ion exchange resin, that permits to obtain demineralized water from tap water.

The ion exchange resin contained in the Demineralizer Bottle is provided with an indicator. When the resin is exhausted, the indicator will change from green to blue. The Demineralizer Bottle must be replaced.

This Demineralizer Bottle may be used to prepare DI water for testing, rinsing or any general use that requires DI water.


INSTRUCTIONS

READ THE ENTIRE INSTRUCTIONS BEFORE USING THE KIT


- Remove the cap and fill the Demineralizer Bottle with tap water.



- Replace the cap and shake gently for at least 5 minutes. The demineralized water is now ready.

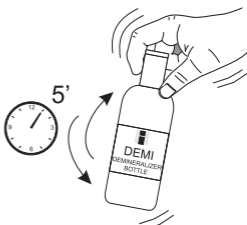

- Flip open the top of the Demineralizer Bottle cap. Pour out the deionized water by gently squeezing the bottle.

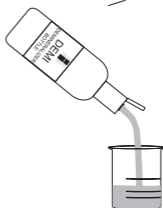

- When the resin is exhausted (its color turns from green to blue) the Demineralizer Bottle must be replaced.

ACCESSORIES

HI 740036

100 mL plastic beaker (6 pcs).

HI 740034

Cap for 100 mL plastic beaker (6 pcs).